


BIOCHROM ANTHOS MULTIREAD400

TECH TIPS: TRANSFERRING DATA TO A PC USING KONTROL+




1. Go to main menu in standalone software on MultiRead 400.
2. Connect instrument to a PC using the serial port or a USB to serial adaptor.
3. Install **Kontrol+** from the CD sent with the instrument.
4. Open **Kontrol+**.
5. Go to **Files>Serial Communication** to establish a connection between the program and the instrument. Set COM port
6. Select **Files>Arange Plates** to download data from the instrument.

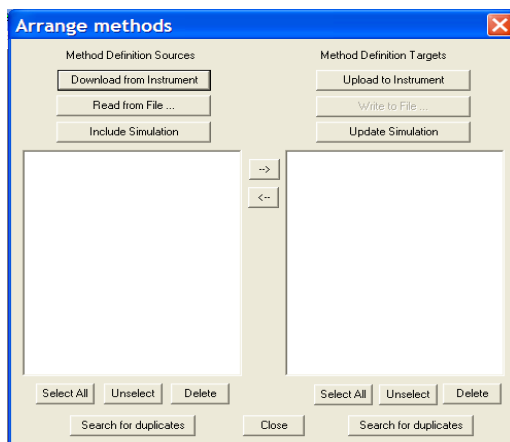


- a. Select **Download from Instrument**.
- b. Select plate or method of interest and select **Write to File**. Save with file name of choice.
- c. Select **Close** to return to the main menu

7. Go to **Special/Printing options....**

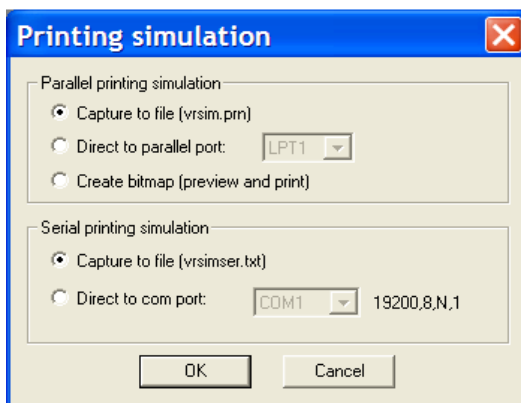

- a. Under **Parallel printing simulation** and **Serial printing simulation** select **Capture to file**.
- b. Select **OK** to return to the **Main Menu**.

9. The features of Kontrol+ mimic the keypad on the instrument. Use the keyboard arrow keys or the mouse on your PC to select **RECALL Plate Data** and confirm by double clicking or pressing enter.

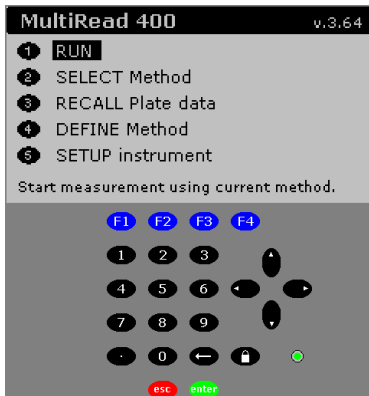

- a. Select plate of interest using the arrow buttons and F4, **Select**.
- b. Select F3, **Print OD**.

10. In analysis program of choice, open **Kontrol+ser.txt** in **Kontrol+** folder in Program Files.

11. When prompted use tab to delimitate data. Now select **Next>Finish**. Data will paste as a matrix into the data analysis software of your choice. Save resulting file.